


## 15ml Centrifuge Adapters for 5ml and 7ml Tubes

The adjustable centrifuge adapters are ideal for properly supporting both 7ml (13 x 100mm) and 5ml (13 x 75mm) tubes in a 15ml centrifuge rotor.


7ml Tube  
(13 x 100mm)



Align the two adapter parts as shown:

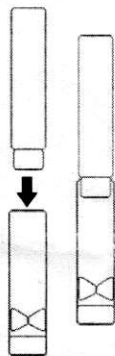

Insert the tube into the adapters as shown:

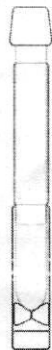


7ml Tube Configuration

5ml Tube  
(13 x 75mm)



Align the two adapter parts as shown:

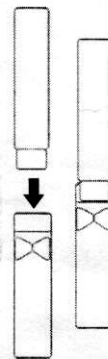

Insert the tube into the adapters as shown:



5ml Tube Configuration