


# Quick Guide

## Monkey NGAL ELISA Kit (KIT 045)

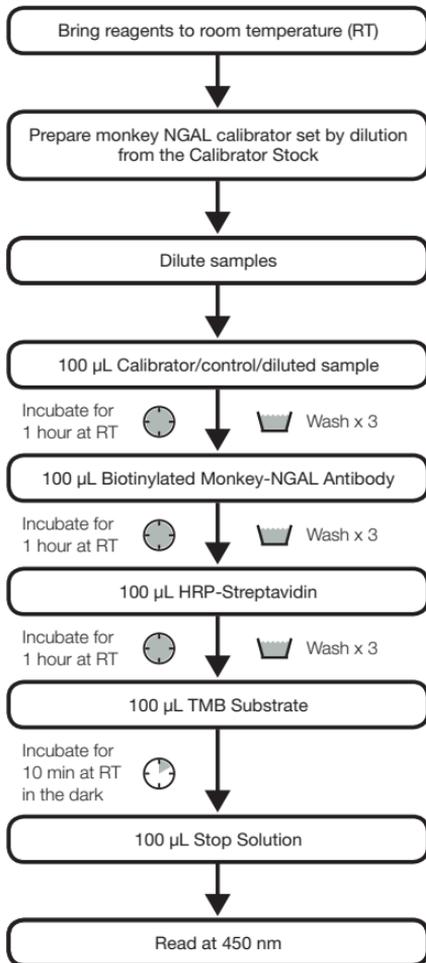


**BIOPORTO**<sup>®</sup>  
Diagnostics


# CALIBRATOR DILUTION GUIDE\*

Monkey NGAL Calibrator Stock

5 mL 1x sample diluent

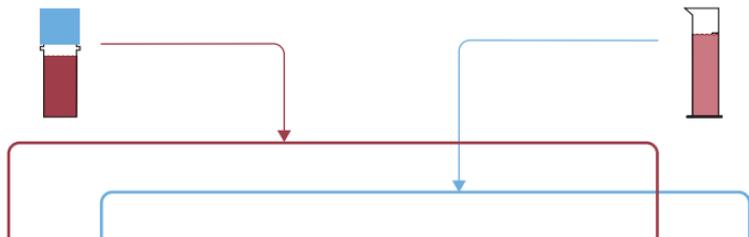

Dilution Factor	1	2	4	10	20	40	100	-
Monkey NGAL	200 $\mu$ L	200 $\mu$ L	100 $\mu$ L	40 $\mu$ L	20 $\mu$ L	10 $\mu$ L	10 $\mu$ L	-
Sample Diluent	-	200 $\mu$ L	300 $\mu$ L	360 $\mu$ L	380 $\mu$ L	390 $\mu$ L	990 $\mu$ L	400 $\mu$ L

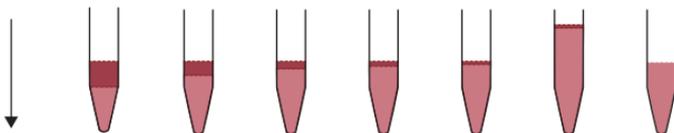

Calibrator #	8	7	6	5	4	3	2	1
NGAL Concentration	1000 pg/mL	500 pg/mL	250 pg/mL	100 pg/mL	50 pg/mL	25 pg/mL	10 pg/mL	0 pg/mL

	↓	↓	↓	↓	↓	↓	↓	↓
Transfer to wells	100 $\mu$ L/ well x 2							

\*Calibrators are to be freshly diluted immediately prior to use.